

Solutions for Your TOUGHEST  
MIXING Applications in

# PHARMACEUTICALS



Assay of Pharmaceutical  
Products

# Assay of Pharmaceutical Products

Assay of products such as pharmaceutical suspensions and tablet coatings is a critical part of the Quality Assurance process. In some cases every batch of each product has to be tested, requiring 100% repeatability of results.

## The Process

Preparation of samples may involve disintegration of solids such as tablets, dispersion of powders into liquids, intimate mixing of liquid/liquid or liquid/solid components to accelerate reactions, etc. Typical process requirements include:


- Efficient particle size reduction and homogenization of the sample before testing is essential in order to obtain 100% consistent and repeatable results.
- The mixing process must be easily and accurately validated.
- Some methods of disintegrating solids, especially on a small scale in the laboratory are difficult to validate; this can also lead to inconsistency between batches.
- Grinding equipment may be difficult to clean, raising potential contamination/hygiene issues.
- The mixer must be capable of performing a wide range of tasks depending on the product being tested, for example disintegration of tablets, dispersion of powders into liquids, etc.
- In some cases a single tablet is tested. The mixing apparatus must be capable of processing such small volumes with the same efficiency as larger samples.

# The Solution

A Silverson High Shear Laboratory Mixer can carry out these duties with great efficiency and repeatable results. Operation is as follows:


## Stage 1

The vessel is charged with appropriate base fluid. The mixer is started and the solid/powdered ingredients are added. The powerful suction created by the high speed rotation of the rotor draws liquid and solids into the workhead where they are rapidly disintegrated.

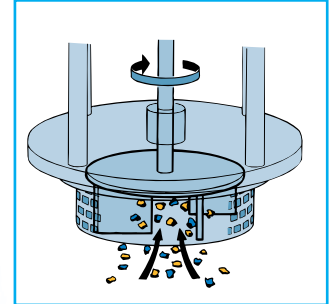


## Stage 2

The materials are subjected to intense but targeted high shear in the precision machined workhead before being forced out through the stator and circulated back into the mix. Simultaneously fresh material is drawn into the workhead.

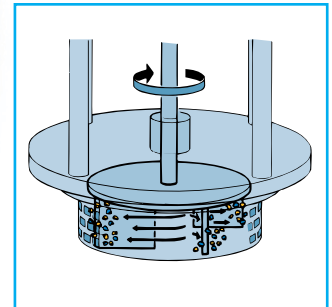

## Stage 3

In a short mixing cycle the entire contents of the vessel passes through the workhead many times, ensuring a completely homogenous mix is obtained.

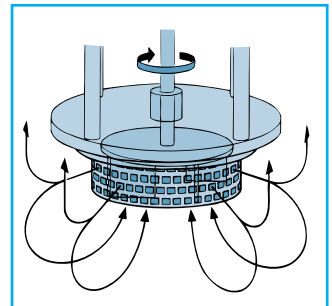

# The Advantages




The Heavy duty L5M-A laboratory mixer has a powerful 1 hp (0.75 kW) motor with variable speed control up to 10,000 rpm. The mixer features a programmable digital timer, ammeter and tachometer, allowing accurate validation of how products perform during processing. This offers a number of advantages:

- 100% repeatable results.
- Easy validation of process.
- Faster mixing reduces risk of degradation of active ingredients.
- Over 40 interchangeable mixing assemblies are available, offering unrivalled versatility.
- Silverson offers the facility to modify mixers to suit client's individual requirements.



In addition to the standard L5M-A model, Silverson offers a number of additional models for this application:

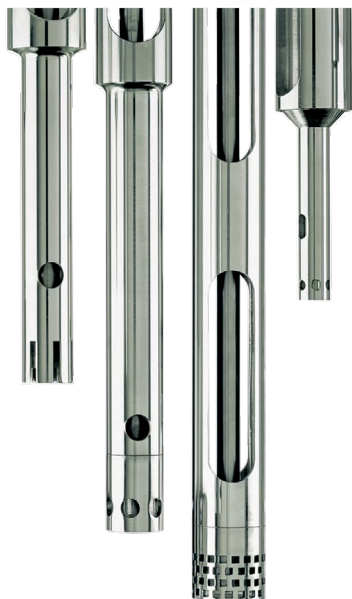

Tubular Mixing Assemblies



Sealed Unit Laboratory Assemblies

- **Tubular Mixing Assemblies**

A series of interchangeable tubular assemblies suitable for use in narrow-necked containers is available for volumes from 1 - 500ml.

- **Sealed Unit Laboratory Mixers**

Designed for processing highly infective or sterile/aseptic substances. Accepts 500ml and 1000ml glass vessels and stainless steel containers up to 5000ml. Flexible tubular assemblies available for small volumes - 225ml, 25ml and 7ml.

- **Further Modifications**

Silverson can offer optional modifications to customise units to the client's requirements. These include:

- A modified bench stand arrangement to ensure that the rise and fall mechanism returns the mixing assembly to the same position in the vessel each time.
- Special rotor/stator assemblies suitable for tablet disintegration can be offered, including assemblies and vessels capable of disintegrating a single tablet.
- Temperature probe with integral digital readout.
- USB port connecting the mixer to a computer for use with the Silverson "Datalogger" program. This records speed and power draw over time and displays it as a graph in real time. The information is also saved in a database.



For more information click here to go to [www.silverson.com](http://www.silverson.com)

Silverson Machines, Inc. 355 Chestnut Street, East Longmeadow, MA 01028  
Ph: (413) 525-4825 [sales@silverson.com](mailto:sales@silverson.com)

Information contained in this report is believed to be correct and is included as a guide only. No warranty is expressed or implied as to fitness for use or freedom from patents. Silverson Machines reserve the right to change product specification without notice.

Issue No. 54PA4